

LABOPROTECT – the protective air exhaust system

LABOKLAV can be equipped with an exhaust air filter. It is suitable to be used in laboratories sterilising pathogen material. The exhaust air filtration system corresponds with the valid regulations. Condensate is being held back in front of the filter, dropping back into the sterilisation chamber and is heated out during the cycle. Integrated into the hydrophobic 0,2 µm filter element is a PT 100 temperature sensor to control that the filter element is completely sterilized during the cycle. Integrated into the housing of the steriliser the filter element is easy to reach and can be replaced within minutes by the user. A stainless steel housing protects the filter element. A filter counter is integrated into the control. The control will remind the user to replace the filter in time.

