

LABOSAFE - protection system for drain pipes

- exhaust steam condenser included into every unit
- condensate temperature adjustable

The protection of the laboratory drain pipes from hot steam is included into the basic version of Laboklav. The discharge system of Laboklav connects steam, condensate and water outlet into one discharge pipe. Tap water is mixed with the condensate leaving the autoclave. By using a PT 100 temperature sensor and the possibility to adjust the condensate and steam drain temperature in the autoclave parameter settings the unit can be adjusted to the type of pipes and the temperature stability of the pipes used in the laboratory.

